

LabSolutions Direct Setup Guide

Introduction

“LabSolutions Direct” enables you to control LC systems (incl. setting parameters, running analysis, and system monitoring) from your smartphone, simply by adding a wireless router to the existing system (*1).

Note that settings are required to enable/disable the LabSolutions Direct feature. Complete the settings by referring to this Setup Guide (*2, *3).

*1 “LabSolutions Direct” does not support controlling/monitoring GC instruments.

*2 Enabling this feature allows the computer to receive external communications. Please only use this feature in an intranet environment where security is guaranteed.

*3 “LabSolutions Direct” supports Windows 7 Professional 32bit/64bit.

Supported Devices


iPhone 4S / iPhone 5 / iPod touch (4G) / iPad (1-4G), and Android 2.3-4.0

* LabSolutions Direct may not operate properly on some Android devices.

Windows is a registered trademark of Microsoft Corporation in the United States and other countries.


iPhone, iPod touch, and iPad are trademarks of Apple Inc.
Android is either a trademark or registered trademark of Google Inc.

1. Enabling the feature

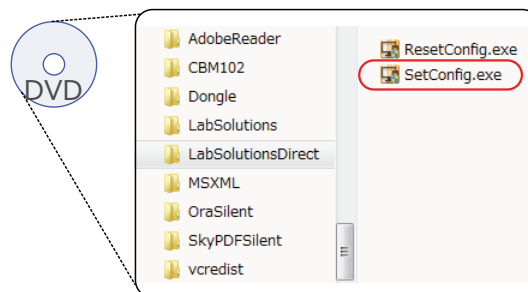

- ① Execute the "SetConfig.exe" included in the installation media.
- ② Click OK in the confirmation box to complete the settings. No more settings are required at the next startup.

2. Disabling the feature

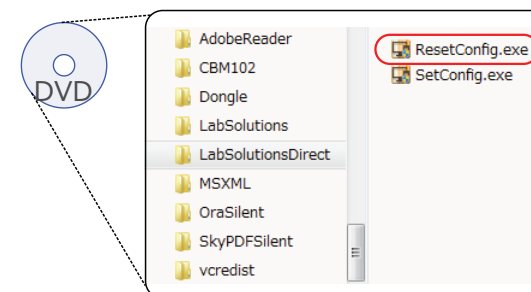

- ① Execute the "ResetConfig.exe" included in the installation media.
- ② Click OK in the confirmation box to complete the settings.

For the operation procedure of LabSolutions Direct, refer to the “LabSolutions Direct Quick Guide,” located at “C:/LabSolutions/Manual/LabSolutionsDirectQuickGuide.pdf” (when LabSolutions is installed on the C drive.)