

Preparation for Demonstrating i-PeakFinder

For the LabSolutions DB/CS, the following preparation is necessary to use the demo data.

1. Create a demonstration project

- (1) On the [Administration] tab of <LabSolutions Main>, Click [Project Administration] to start it.
- (2) Click the [Add] button,
With project name "Demo", create a project.

2. Select the project [Demo]

- (1) On <LabSolutions Main>, select the [Postrun] tab.
- (2) Click the [Select Project] button of <LabSolutions Main>
Select the project name "Demo".

3. Open the browser and copy the demo data

- (1) On the [Postrun] tab of <LabSolutions Main>, [Browser] Click to start.
Then the following folder will be created automatically.
"c:/LabSolutions/Data/Demo"
- (2) Copy all demo data (files under "c:/LabSolutions/Demo/i-PeakFinder")
to the above folder.
* For copying, use Windows Explorer.